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FUNCTION

TOF portable eye wash units with water tank and hand held eye washers comply with CE - DIN 12899 - UNI 9608 standards and they feature a high water flow, but a soft water spout for an effective decontamination of injured body parts, which have accidentally been hit by acid or corrosive chemicals or another harmful substance. Eye washers are indispensable in all areas at risk. They are easy to recognize by their green or red colour and a pictogram sign complying with the standards.

INSTALLATION AND OPERATING INSTRUCTIONS

The tank is equipped with a hand held eye washer / body shower (DIN 12899 - UNI 10271) delivering a soft and foamy water spout. The eye washer is supplied with soft eye protection cups and anti-dust covers.

The valve is opened by slightly pressing the hand lever equipped with a stop device that keeps the valve open and which allows a hands free rinsing process.

TECHNICAL DETAILS FILLING THE TANK

The tank has to be filled with 15 litres of drinking water and pressurized with air at 6 bar.

WATER FLOW

6 l/min for the eye-washer (according to the standards).

WATER TEMPERATURE: 15°C - 35°C

Warning: excessive water temperatures might accelerate some chemical reactions.

In this case we suggest to ask a specialist for advice.

MAINTENANCE

According to the local standards of the country where the eye washer is located.

REPAIRS

Eye washers and body showers may only be repaired in our premises or by skilled and qualified staff. TOF does not undertake any responsibility for alterations or repairs without its consent.

WARRANTY

TOF gives warranty in accordance with the sales contract and the general sales conditions. All products carry a warranty on assembly and manufacturing faults during six months after the delivery date indicated on the shipping documents. The warranty is strictly limited to faults discovered under normal circumstances and correct use. No warranty claim will be accepted for defects caused by alterations or incorrect use by the customer.

IMPORTANT

These instructions can be applied to all TOF eye washers of series 1100.

The drawings do not always show the supplied item.

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